

PRODUCT DATA SHEET

LABTACTICS WASHERS

The Labtactics BOD washer offers features that more expensive units have but for less money. They include; washing, sanitizing, rinse agent, and DI rinse of both BOD & other glassware. Our custom designed spray arm and rack leaves BOD bottles sparkling clean. Sanitizing is accomplished by heating the water to 160 F / 71 C during the rinse cycle, and incorporating DI water for the final rinse. By supplying our washers with two separate glassware racks you are able to wash numerous types of glassware. Our BOD rack allows the operator to quickly & easily insert or remove their BOD bottles from the rack. Another great feature is that our racks can be used to transport your glassware. By lifting out the BOD or lab glassware rack with bottles, you can move them where you like or simply set them to the side for later



use. Install bottle washer on the bench or on rollaway cart, attach quick disconnect fill/drain hose to faucet and turn on hot water. Push DI hose onto DI valve, fill the reservoir with DI water then turn the DI valve on. To start washing just add soap and BOD bottles. In approximately 25 minutes you will have sparkling clean DI rinsed BOD bottles. 115V/60Hz.

LT716
BOD Washer
White Steel Case, Plastic Door Front
22 ½" x 19 ½" x 17 ½"
304 Stainless Steel
16 ½" x 15 ½" x 12"
16 BOD Bottles
Plastic / 6061 Aluminum
16" x 15" x 10"
Rubber Coated Wire / Plastic
115 Volt / 60 Hertz
1.4 Amps
10.4 Amps
160 Degrees F (for 7 minutes)
Hard or Soft Water 50 to 150 Degrees F (soft water is always best)
20 – 90 Psi
Light/BOD 1.6US/gal., Medium 2.4US/gal., Heavy 3.2US/gal.
.8 US/gal.
Yes
Yes for Powder Detergent
90 Days
76 Lbs.
58 Lbs.
Washer 25" x 22" x 21" / Standard Rack & DI Reservoir 18" x 20" x 19"

Labtactics Specifications

Custom washer available, please supply test tube dimensions or part number. Call for quotes

www.garvermfg.com

765-964-5828

sales1@garvermfg.com